



Uniblitz[®] LS2

2mm Uni-Stable Laser Shutter

Overview

The Uniblitz LS2 is a high-performance laser shutter, ideal for precision exposure control, laser switching, and other applications that benefit from the precise, repeatable characteristics of the LS shutter series. These include low level laser chopping, and pulse gating, selection, and modulation to 400Hz. Uni-stable shutter devices, like the LS2, require power to hold the blades in the open state (though the device can be custom-configured for "normally open" operation).

Specifications

Electrical Specifications	
Coil resistance	48 OHMS
Voltage to Open	+65 VDC
Hold Voltage (Nominal) ¹	+10 VDC

Key Features

- 2mm aperture
- Laser Energy Rating: 5 W/mm² (with "ZM" blades)
- Configured for the <u>VCM-D1</u>
 <u>Shutter Driver</u>
- RoHS Compliant
- Transfer time on opening:0.3 milliseconds
- Transfer time on closing:
 0.5 milliseconds

Mechanical SpecificationsWeight Housed210.0 gOperating Temp.0 - 80 °CMax. Opening Bounce15%Max. Closing Bounce5%Max. Freq. of Operation 2100 Hz / 400 HzNumber of Shutter Blades1

¹ Voltage level required across actuator coil when being held in the open position.

² (Continuous/Burst) Continuous frequency rating specified at shutter's minimum exposure pulse. Burst frequency rating specified for four (4) seconds maximum with one (1) minute minimum between bursts.



Need Support? Please visit our website or email us at **info@uniblitz.com**. Tel: +1 585-385-5930 | Fax: 585-385-6004 | 803 Linden Ave. Rochester, NY 14625 Updated 5/24 | Document Version 6.0 | ©2024 Vincent Associates

Uniblitz[®] LS2

Shutter Timing Data



¹ Under no circumstances should any type of lubricant be applied to the shutter blade area. Lubricating the shutter blades will likely slow the shutter down and may eventually render it inoperable.

LS2 (w/ VCM-D1 and "T" blades) ¹ Time (msec.)		
0 - A	Delay time on opening after current a	applied 0.7
A - C	Transfer time on opening	0.3
0 - C	Total opening time	1.0
C - E	Min. dwell time with min. input pulse	0.7
B - F	Min. eqivalent exp. time	1.1
E - G	Transfer time on closing	0.5
A - G	Total window time	1.5
MET	Min. exposure time	1.0
ТЕР	Typical exposure pulse	>1.7

Product Options

LS2 23456-7-8

Ex: LS2S2ZM0-EC-21



² Other blade coating options may be available by special order.

³ Input side only; Teflon[®] coating is on opposite side to protect shutter blade surface. Light source must be input to the reflective side only.

⁴ Mounts are only compatible if #2 housing is optioned as well.

Uniblitz[®] LS2

Electronic Sync.

The electronic synchronization system provides a feedback signal (through the driver utilized) after the shutter transfers to the open state. The system incorporates an infrared emitting diode, an infrared sensitive detecting transistor, and an interrupting vane. The vane is attached to the shutter so as to block the light path between the emitter and detector in the closed position. When the shutter transfers to the 80% open position, the vane is removed from the infrared light path, allowing the emitter to switch the detector to the active state. For the LS2, this system uses a similar activation flag attached to the mechanism, which triggers a reflective emitter/ detector device. No connection to the designated synchronization pins when no electronic sync. is selected.



Uniblitz® LS2 Technical Drawings

Un-Housed



Connector Layout







Uniblitz® LS2 Technical Drawings

501A-S7 Adapter Cable¹ (Not included)



¹ 501A-S7 Shutter Adapter Cable is required for use with VMM-D3, VMM-D4, D880C and any VCM-D1 purchased prior to November 2022.